ULINE

MOBILE FLAMMABLE STORAGE CABINET

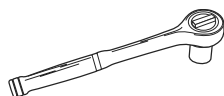
1-800-295-5510
uline.com

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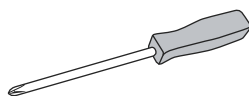
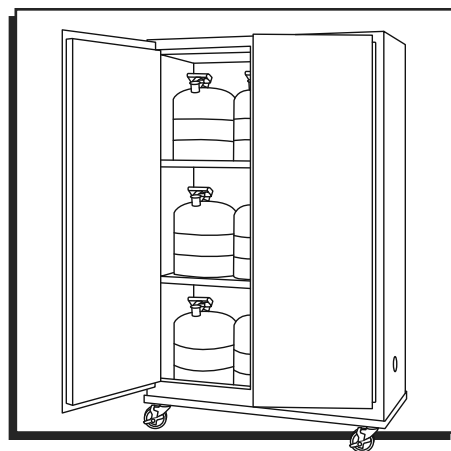
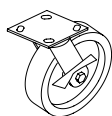
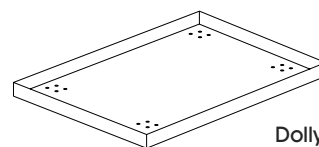
Dolly Tray x 1

CABINET DOLLY ASSEMBLY

1. Align caster plate with bolt holes on dolly tray. Insert four bolts through dolly tray into caster plate. Attach bolts with lock washers and nuts. (See Figure 1)
2. Tighten nuts with a 14 mm wrench.
3. Repeat for remaining casters.


NOTE: Before setting storage cabinet on dolly, ensure all contents are removed from cabinet.

IMPORTANT! Weight capacity of the dolly is 1,000 lbs. This includes weight of storage cabinet and contents.

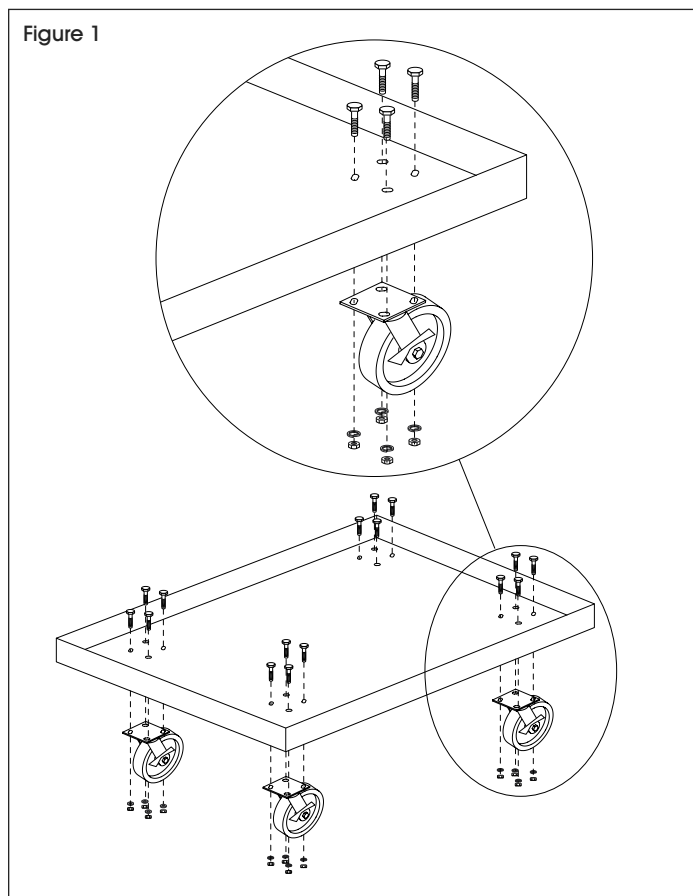Figure 1


CABINET TO BASE ASSEMBLY

NOTE: Before setting storage cabinet on mobile base, ensure all contents are removed from cabinet.

1. Engage brakes on all four casters of mobile base.
2. Carefully lift cabinet onto mobile base, ensure cabinet rests within outer lip of base. (See Figure 2)

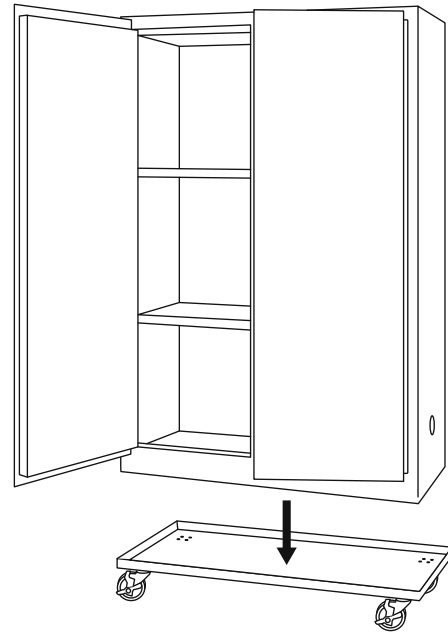Figure 2


SELF-CLOSING CABINET INSTRUCTIONS


1. Install fusible link assembly with screw. (See Figure 3)

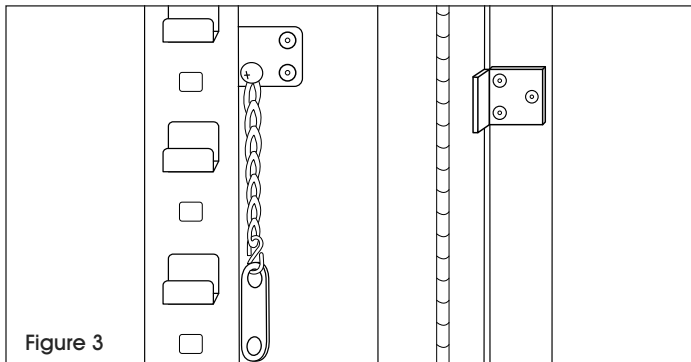NOTE: Brackets are located halfway down inside of right-hand door.

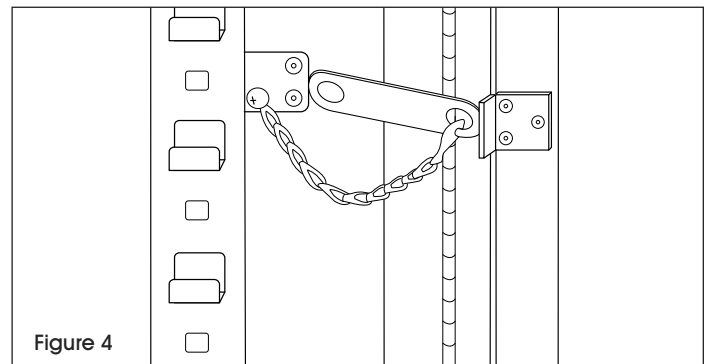2. To hold doors open, prop link between brackets. (See Figure 4)

CABINET GROUNDING INSTRUCTIONS

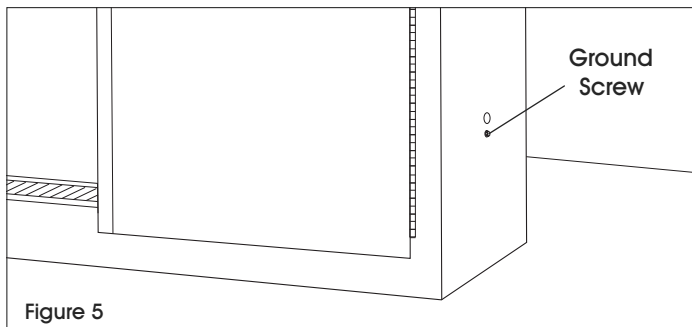
NOTE: To ensure safety in facilities, it's important to eliminate the chance of an electrostatic discharge while handling flammable fluids. Refer to NFPA 77 for further information.

CAUTION! If dispensing Class 1 flammable liquids from a container held within safety cabinet, ground the cabinet. It is recommended to ground the cabinet when dispensing Class 2 combustible liquids if the liquids are near, at or above the liquid's flashpoint temperature. Be aware of these external and internal grounding considerations.

EXTERIOR CABINET GROUNDING


1. Locate the safety cabinet's ground screw just below the green dot on the right-hand side. (See Figure 5)


2. Use a 5/16" (8 mm) wrench or nut driver to loosen screw. Attach the ground wire and retighten screw. (See Figure 6)

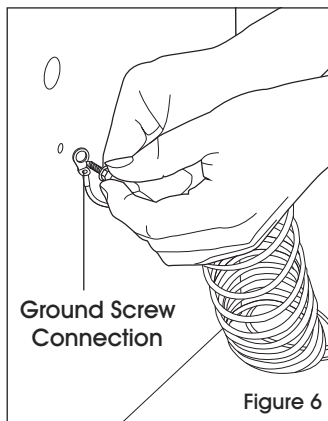NOTE: Ground wire not included.

3. Attach opposite end of ground wire to a ground bus or ground rod. Ground cabinet to earth. (See Figure 7)

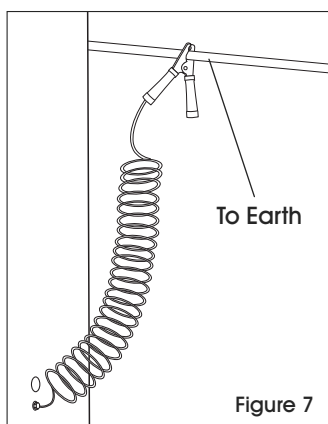
INTERIOR CABINET GROUNDING

When dispensing liquids inside a safety cabinet, it is important to ensure good electrical continuity through metal-to-metal contact of ground wires.

1. Bond between safety can and drum or between drum and upper vent hole. (See Figure 8)

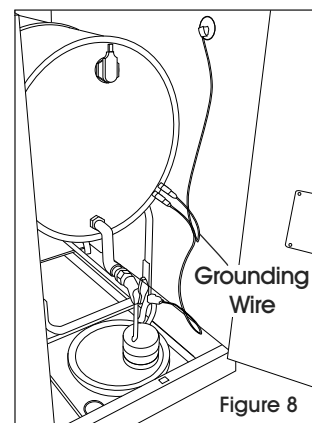NOTE: User may have to remove some paint to contact metal. If necessary, sand off a small area of paint where clamp is attached to ensure clamps contact metal.

RECOMMENDATIONS:

- Use commercially available grounding wires, as they are made for industrial service duty.
- Ensure all connections are secure.
- After making connections, user should discharge by touching ground before opening containers.
- Follow these guidelines to help comply with the law when dispensing Class 1 flammable liquids, and help eliminate the chance of a static electrical discharge and the dangerous ignition of vapors.

NOTE: This information was prepared by the manufacturer and is provided as a service to our customers. Uline is not responsible for the accuracy or completeness of this information.

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