

ULINE H-5100 FAUCET MOUNT EYEWASH STATION

1-800-295-5510
uline.com

TOOL NEEDED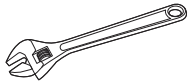

Adjustable Wrench

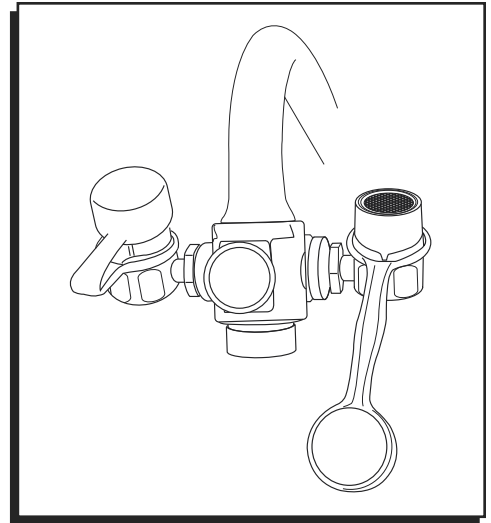INSTRUCTIONS

INSTALLATION

1. Remove present aerator or other fitting from faucet outlet.
2. Mount the eyewash station to faucet spout. The eyewash station inlet has a 55/64"-27 female thread.

NOTE: Due to the variety of faucets, it may be necessary to use an adapter to mount the eyewash station on the faucet spout. Three adapters come with the unit: a larger adapter with 15/16"-27 thread size to fit Delta and American Standard faucets, a smaller adapter with 13/16"-27 thread size to fit Kohler and other faucets, and a 3/8" IPS adapter to fit Water Saver faucets. An additional 13/16"-24 female adapter is available that fits Chicago brand faucets.

IMPORTANT! When mounting eyewash station to faucet, be careful not to overtighten adapter.

3. If necessary, use a 13/16" or adjustable wrench to adjust the angle of the outlet heads to obtain the desired angle for the eyewash water flow.
4. Attach the enclosed sign to wall or backsplash behind sink.

OPERATION

1. Turn on cold water side of faucet.

IMPORTANT! Use only cold water when eyewash station is in operation. Hot water may cause scalding or other eye injury.

2. Pull knob on eyewash station unit forward to activate the eye wash. Water pressure will hold the eyewash in operation, leaving the user's hands free.
3. To return to normal faucet operation, push knob back to original position. When the faucet is turned off, the knob will return to original position automatically.

MAINTENANCE

- Periodic cleaning of the eyewash aerators is recommended to ensure proper water flow.
- Keep plastic covers on outlet heads when the unit is not in use.
- The eyewash station should be tested weekly.

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