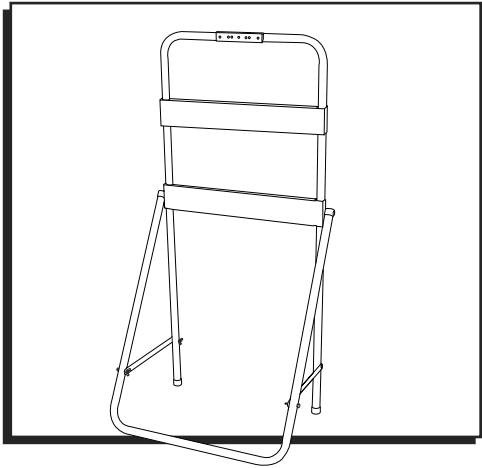
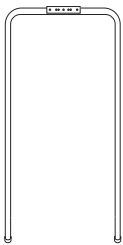
ULINE H-5193 EYEWASH STAND

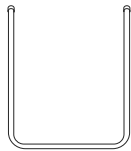
1-800-295-5510
uline.com


PARTS


Back Leg x 1

Front Leg x 1

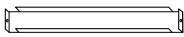
Back Brace x 2

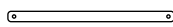Leg Brace x 2

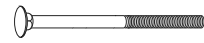1 3/4" Cross Head Bolt x 9

3/4" Carriage Bolt x 2

Washer x 11

Lock Nut x 4

Wing Nut x 7

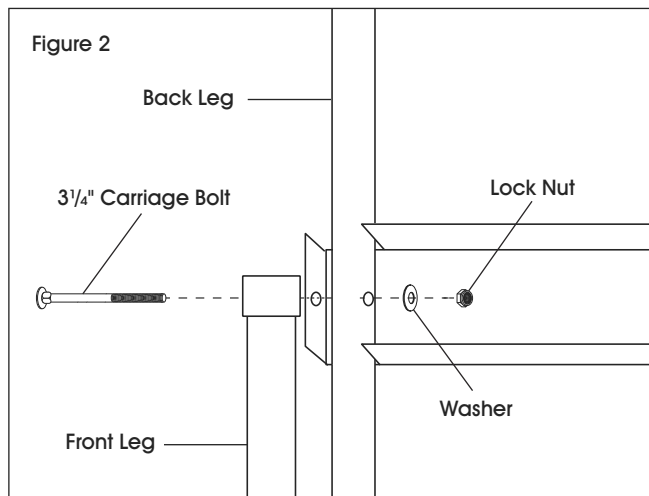
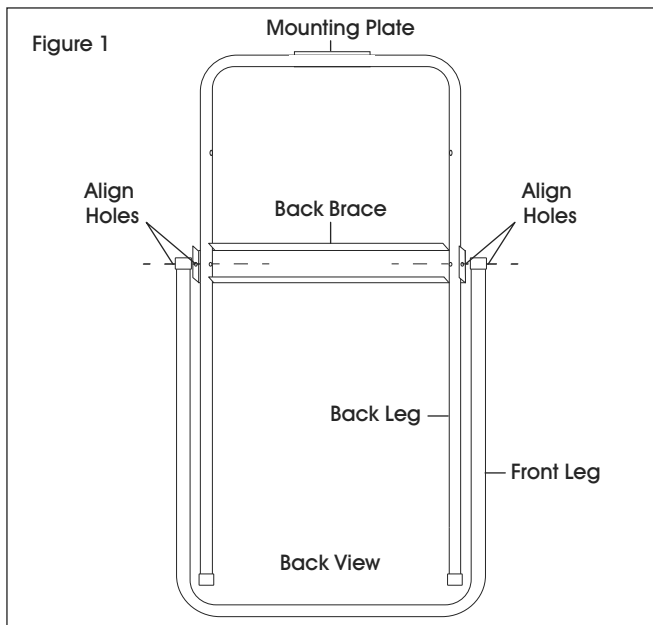
ASSEMBLY

1. Align middle holes of the back leg with holes of the back brace. (See Figure 1)

 **NOTE:** Flat panel of the back brace should face the same direction as mounting plate.

2. Align the holes of the front leg with holes of the back brace and back leg. (See Figure 1)


3. Secure the back brace, back leg and front leg by using a 3/4" carriage bolt, washer and lock nut. Repeat for opposite side. Do not over tighten. (See Figure 2)


 **NOTE:** Bolts enter through the outside of the unit and are fastened from the inside.

ASSEMBLY CONTINUED

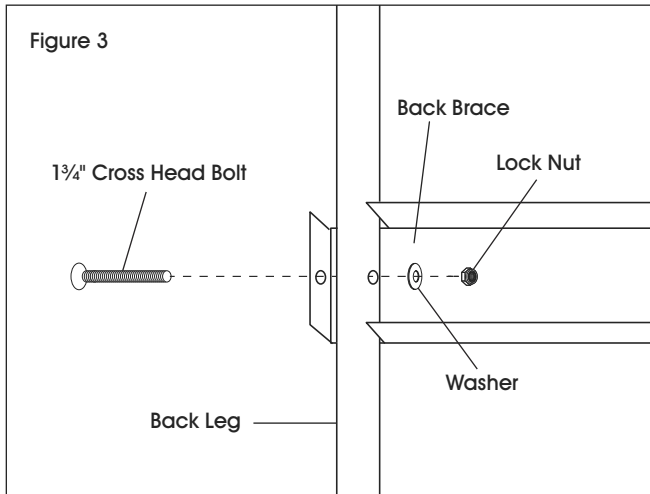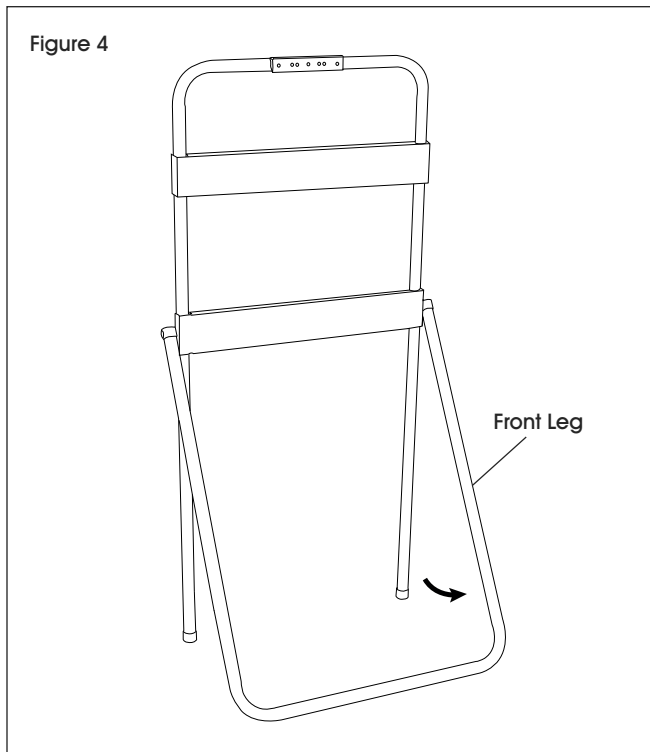
- Align side holes of remaining back brace with top holes of the back leg.
- Secure back brace to the back leg by using a 1 $\frac{3}{4}$ " cross head bolt, washer and lock nut. Repeat for opposite side. (See Figure 3)

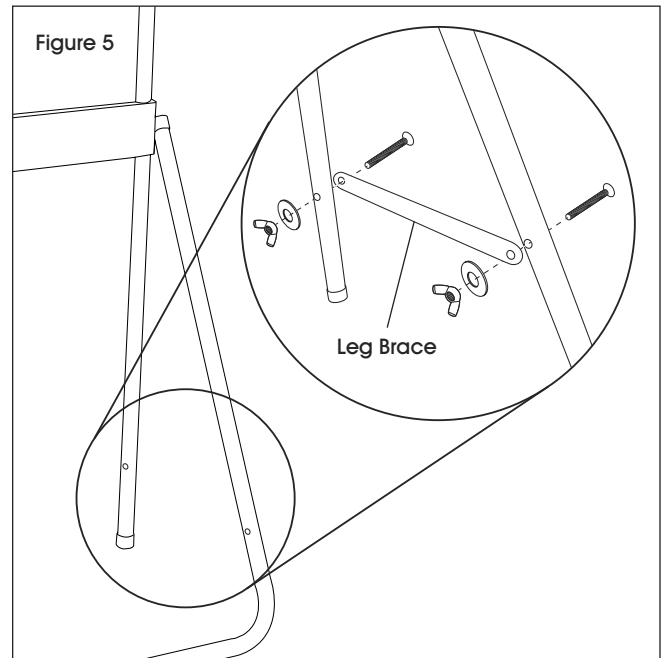- Pull the front leg out from the unit as shown. (See Figure 4)

- Attach one leg brace to back leg and front leg using two 1 $\frac{3}{4}$ " cross head bolts, two washers and two wing nuts. Repeat for remaining leg brace. (See Figure 5)

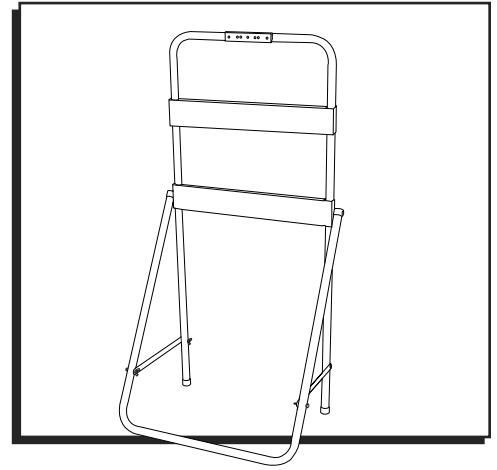
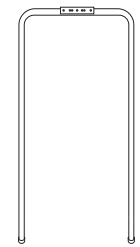
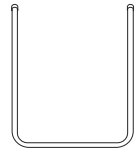
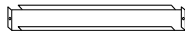
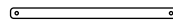
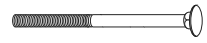
-  **NOTE:** Additional hardware can be used to install eyewash station. Mounting bracket is included with the eyewash station.
-  **NOTE:** When attaching an eyewash station, refer to the installation instructions included with the unit.

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