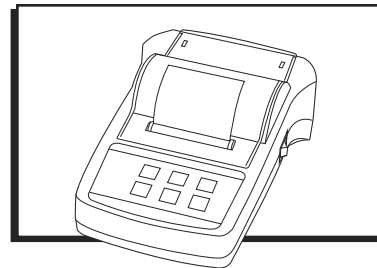
ULINE H-8093

OHAUS SF40A[®]

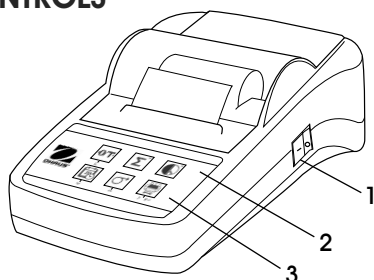
SCALE PRINTER

1-800-295-5510
 uline.com


OVERVIEW OF CONTROLS

CONTROLS


CONNECTIONS

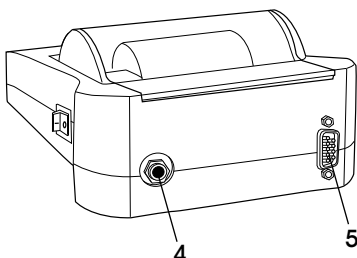CONTROLS AND CONNECTIONS

#	DESCRIPTION
1	On/Off Switch
2	Pilot Lamp – Status Indicator
3	Control Panel
4	Power Connection
5	Interface Connector

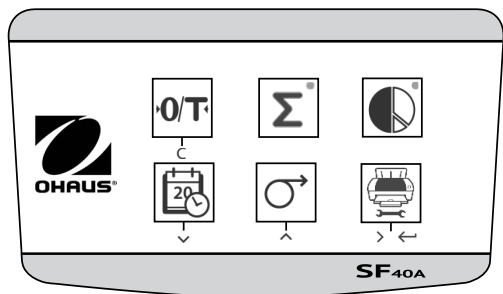CONTROL PANEL

CONTROL FUNCTIONS

KEY	DESCRIPTION	KEY	DESCRIPTION
	Zeroing/Taring the balance Abort – Exit		Paper feed Scroll up menu option
	Printout of date and time Scroll down menu option		Statistics application
	Totalizing application		Printing of stable weight values Opening the menu – select next menu option Save settings

SETUP

SAFETY

The printer has been tested for the connections and intended purposes documented in these operating instructions. It is your responsibility to perform the necessary tests related to the methods and purposes of its intended use. Observe the following safety measures.

CAUTION! The printer must only be used indoors and never in hazardous areas.

- The electronics of the printer are protected against splashing water and dust. However, since the printer housing is not watertight, it should not be used in the presence of liquids.

- Follow cleaning instructions on page 8. Protect power plug against wet conditions.
- Never open the printer housing. If you have any problems with your printer, contact Uline Customer Service at 1-800-295-5510.
- If the equipment is used in a manner not specified by the manufacturer, the protection provided by the equipment may be impaired.

POWER SUPPLY

The printer is supplied with an AC adapter with a country-specific power cable. The power supply is suitable for all line voltages in the range: 100 – 240V AC, 50 – 60 Hz. (See Page 11 for Specifications)

SETUP CONTINUED

CAUTION! First, check the local line voltage is in the range 100 – 240V AC, 50 – 60 Hz and determine whether the power plug fits your local power supply connection. If this is not the case, do not connect the printer or the AC adapter to the power supply. Contact Uline Customer Service at 1-800-295-5510.

- Only plug adapter into a grounded socket.

WARNING! Before operating, check all cables for damage.

- Guide the cables so that they are not damaged or interfere during the weighing process.
- Make sure that the AC adapter cannot come into contact with liquids.
- The power plug must always be accessible.

PRINTER SETUP

- Default printer settings are shown below. (See Figure 1)

Figure 1

Header:	No Header	Baud rate:	9600
Date Format:	DD MMM YYYY	Data bits:	8
Time Format:	24:MM:SS	Parity:	None
Char Set:	CP1252 Window Latin1	Stop bits:	1
Balance Feature:	Balance Feature Enable	Flow control:	Xon/Xoff
O/T Settings:	Perform Zero		

There are two ways to change the default setting:

- Use accompanying software.
- Operate on the printer only.

PRINTER SETUP USING ACCOMPANYING SOFTWARE

1. Attach the interconnecting cable to the printer interface on the computer and to the printer.
2. Connect printer to the main power supply.
3. Use the main switch <I/O> to power on.
4. Insert CD into the computer and install the printer software.
5. Run the software and set up the RS-232 parameter of computer, then click the Connect button. (See Figure 2)

Figure 2

COM Settings			
Please config the COM port setting.			
Port:	COM1	Baud rate:	9600
Data bits:	8	Parity:	None
Stop bits:	1	Flow control:	Xon/Xoff
<input type="button" value="Connect"/>		<input type="button" value="Disconnect"/>	

6. The following window will appear if connection is successful. Click OK to continue. (See Figure 3)

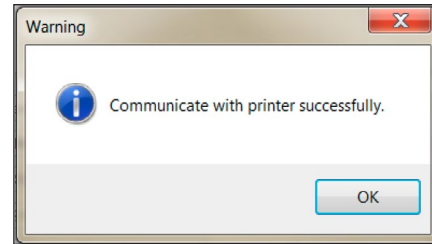
Figure 3

7. The printer Menu setting is now activated and all settings can be changed. (See Figure 4)

Figure 4

Menu Setting			
Date Time:	<input checked="" type="radio"/> Keep Printer Time <input type="radio"/> Current System Time <input type="radio"/> Custom	2012/08/01 09:51:35	
Header:	No Header	Baud rate:	9600
Date Format:	DD MMM YYYY	Data bits:	8
Time Format:	24:MM:SS	Parity:	None
Char Set:	CP1252 Window Latin1	Stop bits:	1
Balance Feature:	Balance Feature Enable	Flow control:	Xon/Xoff
O/T Settings:	Perform Zero		
Description: CP1252 covers Western European languages such as English, German, French, Spanish, Italian and other languages.			
<input type="button" value="Factory Reset"/>		<input type="button" value="Get Status"/> <input type="button" value="Save Status"/>	

a. **Date Time:** Sets date and time for the printer.

- **Keep Printer Time:** Keep existing date and time in the printer.
- **Current System Time:** Get date and time from your computer.
- **Custom:** Set date and time manually.

b. **Header:** Sets the header of the printing content.

- **No Header:** No header for printing content.
- **Date/Time:** Header with date and time.
- **Date/Time/Balance:** Header with date, time and S/N.

c. **Date Format:** Sets date format.

d. **Time Format:** Sets time format.

e. **Char Set:** Sets code page for different languages.

- **CP 1252 Windows Latin1:** Covers Western European languages such as English, German, French, Spanish and Italian.
- **CP 1251 Windows Cyrillic:** Covers languages that use the Cyrillic alphabet such as Russian, Bulgarian and Serbian Cyrillic.

SETUP CONTINUED

- CP437 MS-DOS Latin US: Covers Western European languages such as English, German, French, Spanish and Italian.
- CP852 MS-DOS Latin2: Covers Central European languages that use Latin script such as Polish, Bosnian, Czech, Hungarian, Croatian, Romanian, Serbian and Slovak.
- CP860 MS-DOS Portuguese: Covers Portuguese language.

f. Balance Features:

1. **Enable/Disable a function:** O/T, Statistics, Totalization.
2. **Printer connection status:** LED light will either blink or stay lit.
 - Balance Feature Enabled: O/T, Statistics and Totalization functions are enabled. The printer connection status LED light will blink if the connection with balance is lost.
 - Balance Feature Disable: O/T, Statistics, Totalization functions are disabled. The printer connection status LED will be lit even if the connection with balance is lost.

g. O/T Setting: Set up the function of the O/T button.

- Perform Zero: Perform Zero function when pressing the button.
- Perform Tare: Perform Tare function when pressing the button.

h. Baud rate, Data bits, Parity, Stop bits, Flow Control: Set the parameters of RS-232 setting.

- Settings should be the same as the scale to be connected. For the scale RS-232 settings, please see scale manual.

Factory Reset

Reset printer to factory default setting.

Get Status

Get current setting from the printer.

Save Status

Save the setting to the printer.

8. Click the Save Status button to save the setting and click Exit to finish the setup.

PRINTER SETUP USING PRINTER ONLY

1. Connect the printer to the main power supply.
2. Use the main switch <I/O> to power on.
3. Press and hold the  button to enter the menu list, the printer will print "-----Menu-----".

- Press  to choose which item to change.

- Press  to change the parameter.

- Press and hold  to save and exit.

- Press  to exit without saving.

The printer will print the content to show the result of every operation. See Operation for total menu structure. (See pages 5-8)

INSTALLING THE PRINTER

CONNECTING THE PRINTER TO THE SCALE


Before connecting, make sure the scale and printer have the same RS-232 parameter setting. For scale Baud rate settings, see Recommended Settings on pages 6-7.

1. Attach interconnecting cable to the printer interface of the scale and to the printer.
2. Switch the scale on.

SWITCHING THE PRINTER ON

1. Connect printer to the main power.
2. Use main switch <I/O> to power on. (See Figure 5)

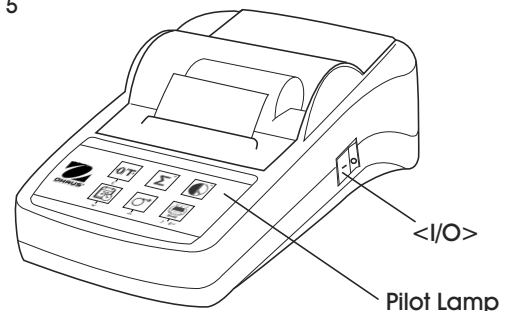Figure 5


3. RS-232 interface only: The pilot lamp will flash until scale is connected. If pilot lamp continues to flash, see Troubleshooting on pages 9-10.

SETUP CONTINUED

INSERTING PAPER

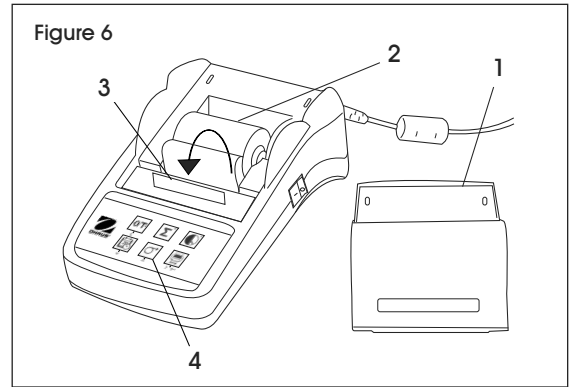1. Pull back of paper cover (1) to remove. (See Figure 6)
2. Insert paper rotary axis through the core of the paper roll (2).
3. Feed paper through the slot in the printing unit in the direction of the arrow (3).
4. Press and hold until enough paper has been fed (4).
5. Replace paper cover (1).


PRINTER TEST

The printer is equipped with a self-test that automatically prints out the character set (standard IBM/DOS).

- Starting the Test – Press the  key and power on the printer. The printer prints out the current settings.
- Terminating the Test – Power off printer.

SETUP MENU

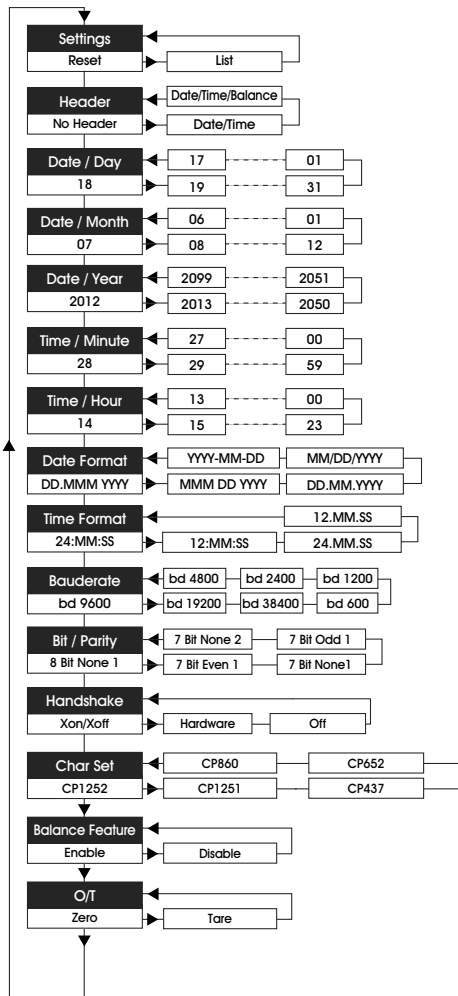SETUP MENU

SAMPLE PRINTOUT




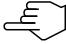
```

-----MENU-----
Settings:
* List

Header:
* No Header

Date Day:
31

Date Month:
07

Date Year:
2012

Time Minute:
00

Time Hour:
18

Date Format:
* YYYY-MM-DD

Time Format:
* 24:MM:SS

Baudrate:
* 9600

Bit/Parity:
* 8 Bit None 1

Handshake:
* Xon/Xoff

Char Set:
* CP1252 Windows Latin 1

Balance Feature:
* Enable





->O/T<- Setting:
* Zero
    
```

* Factory Settings


OPERATION

KEYSTROKE FUNCTIONS

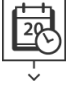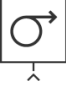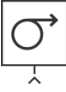OPERATIONAL LEVEL

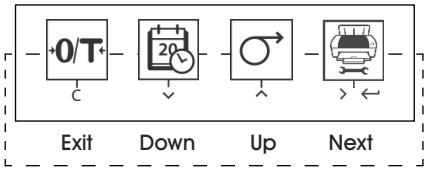 PRESS BRIEFLY	 PRESS AND HOLD
 Zero or tare the balance	
 Print of date and time	
 One line feed	 Continuous feed (until release)
 Print	 Open menu

MENU LEVEL

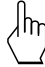
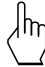
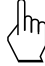
 Exit Down Up Next	 Store
---	--

FUNCTION LEVEL

 Start totaling application	 Conclude totaling application
 Start statistics application	 Conclude statistics application

OPERATION CONTINUED

APPLICATIONS

 PRESS BRIEFLY		 PRESS AND HOLD	
STATISTICS APPLICATION		TOTALING APPLICATION	
 	Start Statistics Application (LED on)	 	Start Totaling Application (LED on)
 	Add a new sample and press 	 	Add a new sample and press 
 	If needed, print subtotal. Press 	 	If needed, print subtotal. Press 
 	Print and end Statistics Application (LED off)	 	Print and end Totaling Application (LED off)

 **NOTE:** Unit change during the application is not permitted.

- Maximum 999 samples.
- Sample weight must be in the range of 70 – 130% of the current average value.
- Minimum sample weights should be larger than 1d.
- During sample change, the scale reading needs to change by more than 100d or the printer will not register the sample change.
- Statistics applications include: number of samples, average value, standard deviation, min., max. and min-max differential and total weight.

RECOMMENDED SETTINGS

CONNECTED DEVICE	DEVICE SETTINGS	PRINTER SETTINGS
DV, AV, PA, AR, SP, TA, NV, NVL, NVT, MB35/45, MB23/25, (T31P, T51P, T71P, T32XW, T32WE, T51XW, T71XW, CKW)*	Printer Mode Baud Rate: 9600 Bit/Parity: 8 bit None 1 Handshake: Xon/Xoff	Factory Settings

OPERATION CONTINUED

RECOMMENDED SETTINGS CONTINUED

CONNECTED DEVICE	DEVICE SETTINGS	PRINTER SETTINGS
EX	Printer Mode Baud Rate: 9600 Bit/Parity: 8 bit None 1 Handshake: Xon/Xoff	Factory Settings
Russian	Printer Mode Baud Rate: 9600 Bit/Parity: 8 bit None 1 Handshake: Xon/Xoff Language: ANSI/WIN Russian	CP1251 Windows Cyrillic
Polish	Printer Mode Baud Rate: 9600 Bit/Parity: 8 bit None 1 Handshake: Xon/Xoff Language: ANSI/WIN Polish	CP852 MS-DOS Latin2
Czech	Printer Mode Baud Rate: 9600 Bit/Parity: 8 bit None 1 Handshake: Xon/Xoff Language: ANSI/WIN Czech	CP852 MS-DOS Latin2
Hungarian	Printer Mode Baud Rate: 9600 Bit/Parity: 8 bit None 1 Handshake: Xon/Xoff Language: ANSI/WIN Hungarian	CP852 MS-DOS Latin2
Portuguese	Printer Mode Baud Rate: 9600 Bit/Parity: 8 bit None 1 Handshake: Xon/Xoff Language: ANSI/WIN Portuguese	CP860 MS-DOS Portuguese
English, French, German, Italian, Spanish, Chinese, Japanese, Korean	Printer Mode Baud Rate: 9600 Bit/Parity: 8 bit None 1 Handshake: Xon/Xoff	Factory Settings

* The Statistic and Totalization functions of printer are supported in the following models: DV, EX, AV, PA, NV, MB23/25. Only the printing function is supported in all models.


OPERATION CONTINUED

RS-232C INTERFACE

The SF40A printer is equipped with an RS-232C interface to connect Ohaus scales. (See Figure 7)

- 9-pin male connector
- Matching to other devices (transmission parameters). (See Connecting the Printer to the Scale on page 3)

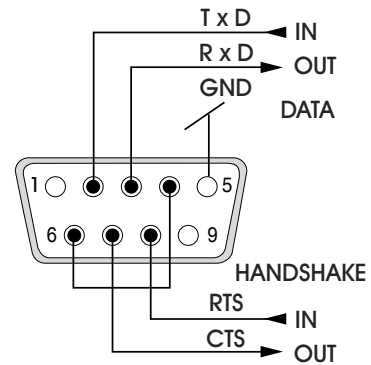Figure 7

MAINTENANCE

Under normal conditions, the printer requires practically no maintenance. If servicing is required, contact Uline Customer Service at 1-800-295-5510

CHANGING THE PAPER ROLL


Insertion of a new roll of paper is described in Inserting Paper on page 4. (For paper roll order number, see Consumables and Spare Parts on page 10)

 **NOTE:** You may obtain the paper rotary axis from the old paper roll.

REPLACING THE RIBBON


1. Remove the paper cover (1). (See Figure 8)
2. Pull paper out of printing unit (2)

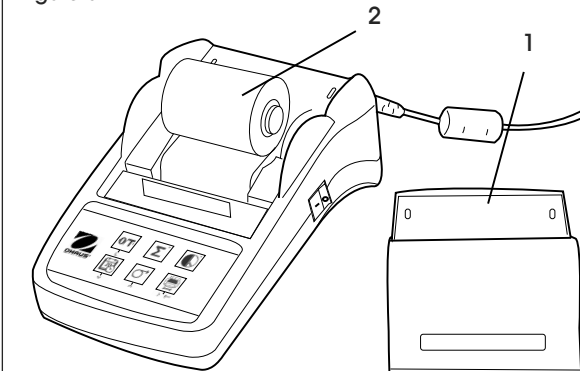Figure 8

3. To remove the ribbon cartridge, press in direction of arrow (3). (See Figure 9)

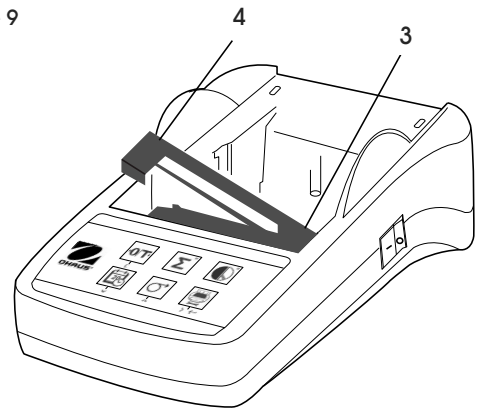Figure 9

4. Insert new ribbon, and if necessary, take up slack with tensioning wheel (4).

 **NOTE:** For the ribbon cartridge order number, see Consumables and Spare Parts on page 10.

 **NOTE:** For insertion of a new roll of paper, see Inserting Paper on page 4.

CLEANING

- Since the printer housing is made of top grade materials, all commercially available mild cleaning agents may be used. (See Safety Precautions on pages 1-2)

TROUBLESHOOTING


OPERATING ISSUE	CAUSES	RECOMMENDATIONS
Printout unreadable.	Ribbon worn or tangled. Service life of printing unit has ended.	Change ribbon cartridge or tense the ribbon properly. Contact Uline Customer Service at 1-800-295-5510.
Pilot lamp does not light.	No power. Power supply fault.	Switch printer on. Contact Uline Customer Service at 1-800-295-5510.
Pilot lamp flashes.	No connection printer to device. Printer unable to recognize device settings for the connected scale.	Make sure that the printer and the scale are connected. Switch the printer off/on. Alternatively change settings of the printer or the connected scale. (See page 4)
----- Timeout -----	Data transfer not completed within 45 seconds.	Move scale to a more stable location. Check the scale.
----- Over Load -----	Balance in overload range. Zeroing out of the zero range.	Check the sample weight. Change the zero range of the balance. Change the setting of the printer from zero to tare.
----- Under Load -----	Scale in underload range. Weighing pan is not in position.	Check the scale.
---- Transmission Error----	Scale does not recognize printer instruction.	Check the interface parameter settings on the printer and scale.
----- Not Executable -----	Scale cannot execute printer instruction.	Check the scale.
--- Date/Time Not Set --- ----- Memory Error -----	Printer error messages.	Check the scale. Reset the printer. Set date/time. Backup battery depleted. Contact Uline Customer Service at 1-800-295-5510.

TROUBLESHOOTING CONTINUED

OPERATING ISSUE	CAUSES	RECOMMENDATIONS
----- Memory Full -----	Statistics or Totaling Memory is full.	Start a new Statistics or Totaling application.
-- Load/Unload Weight --	No sample or the same sample is on the weighing pan.	Put a new sample on the weighing pan or remove it and put sample on the weighing pan again.
Unit change not allowed!	Changed unit during Statistics or Totaling application.	Reset the scale. Change the unit back to the original unit setting.
----- Out of range -----	Sample weight is out of range.	Sample weight must be in the range of 70-130% of the current average value.

If the troubleshooting section does not solve your problem, contact Uline Customer Service at 1-800-295-5510.

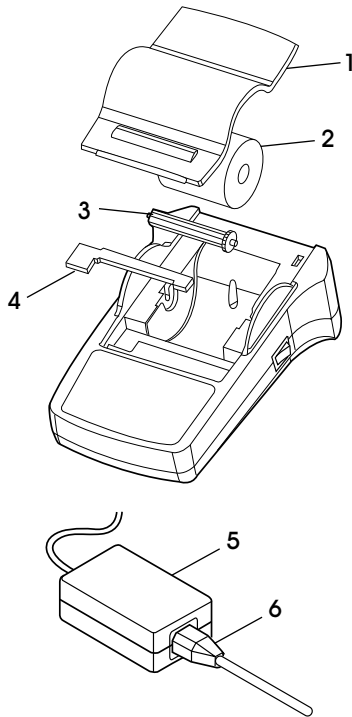CONSUMABLES AND SPARE PARTS

#	DESCRIPTION	QTY.	ULINE PART NO.	MFG. PART NO.
1	Paper Cover	1	-----	30047939
2	Paper Roll	1	S-23229	12120799
3	Paper Rotary Axis	1	-----	30063920
4	Ribbon Cartridge, Black	1	S-23230	12120798
5	AC/DC Adapter (without power cable) 100-240V AC, 50-60 Hz, 12V DC 2.5 A	1	-----	11107909
6	Power Cable CH	1	-----	30064089
	Power Cable EU		-----	12120761
	Power Cable US		H-8093-PCBL	12120762
	Power Cable IT		-----	30064200
	Power Cable GB		-----	12120312
	Power Cable AU		-----	12120313
	Power Cable BR		-----	83033772
	Power Cable JP		-----	12122638
-	RS9 M/F Cable, 1.5 m	1	-----	12122603

SPECIFICATIONS

Power Supply	AC/DC adapter: 12V DC, 2.5 A Primary: 100 – 240V AC, -50 -60 Hz Voltage Fluctuations: Up to $\pm 10\%$ of the nominal voltage Secondary: 12V DC, 2.5 A (with electric overload protection)
Cable to AC/DC Adapter	3-core with country-specific plug.
Power Line Voltage	12V DC
Power Consumption	Max. 2.5 A
Altitude	Up to 2000 m
Ambient Temperature	0° C – 40° C
Relative Humidity	Max. 80% at 31° C, decreasing linearly to 50% at 40° C, noncondensing
Overvoltage Category	Class II
Pollution Degree	2
Printing Unit	Dot matrix printer 5 x 7, 40 character line length
Printing Speed	1.2 lines per second
Ribbon Cartridge	Exchangeable, black
Paper Roll	Standard paper 57.5 x 50-60 mm, integrated in housing, commercial size
Print Quality	Light resistant and thermally stable printing (GLP, GMP, ISO9001)
Interface	RS-232C
Printer Dimensions	(W x D x H) 120 x 203 x 73 mm
Packaging Dimensions	(W x D x H) 255 x 205 x 150 mm
Net Weight	740 g (paper roll included)
Compatibility Limits	Function Date/Time: Not applicable in balances with built-in clock (built-in clock has priority). Functions Zeroing/Taring, Statistics and Totalization: Not applicable on all models of balances.

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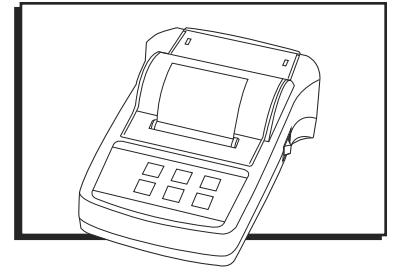
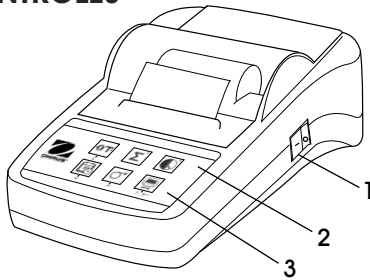
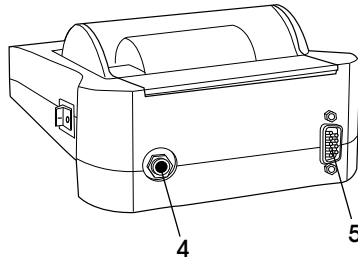
Diagrama 2

COM Settings			
Please config the COM port setting.			
Port:	COM1	Baud rate:	9600
Data bits:	8	Parity:	None
Stop bits:	1	Flow control:	Xon/Xoff
Connect		Disconnect	

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Diagrama 4

Menu Setting			
Date Time:	<input checked="" type="radio"/> Keep Printer Time <input type="radio"/> Current System Time <input type="radio"/> Custom	2012/08/01 09:51:35	
Header:	No Header	Baud rate:	9600
Date Format:	DD MMM YYYY	Data bits:	8
Time Format:	24:MM:SS	Parity:	None
Char Set:	CP1252 Window Latin1	Stop bits:	1
Balance Feature:	Balance Feature Enable	Flow control:	Xon/Xoff
O/T Settings:	Perform Zero		
Description: CP1252 covers Western European languages such as English, German, French, Spanish, Italian and other languages.			
Factory Reset		Get Status Save Status	

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Figure 2

COM Settings			
Please config the COM port setting.			
Port:	COM1	Baud rate:	9600
Date bits:	8	Parity:	None
Stop bits:	1	Flow control:	Xon/Xoff
<input type="button" value="Connect"/>		<input type="button" value="Disconnect"/>	

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Figure 4

Menu Setting			
Date Time:	<input checked="" type="radio"/> Keep Printer Time <input type="radio"/> Current System Time <input type="radio"/> Custom	2012/08/01 09:51:35	
Header:	No Header	Baud rate:	9600
Date Format:	DD.MMM.YYYY	Date bits:	8
Time Format:	24.MM.SS	Parity:	None
Char Set:	CP1252 Window Latin1	Stop bits:	1
Balance Feature:	Balance Feature Enable	Flow control:	Xon/Xoff
O/T Settings:	Perform Zero		
Description: CP1252 covers Western European languages such as English, German, French, Spanish, Italian and other languages.			
<input type="button" value="Factory Reset"/>		<input type="button" value="Get Status"/> <input type="button" value="Save Status"/>	

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