
# ULINE H-9460 EYEWASH DRENCH HOSE – WALL MOUNT

1-800-295-5510  
uline.com


## TOOLS NEEDED

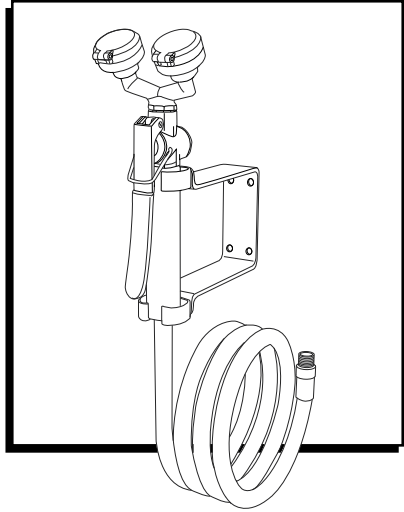Electric Drill


1/4" Drill Bit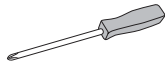

Phillips Screwdriver


## PARTS

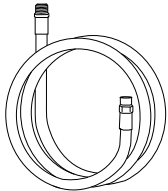8' Hose x 1


Handle x 1

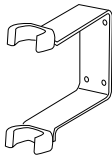
Wall Bracket x 1

Screw x 4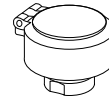

Spray Head x 2

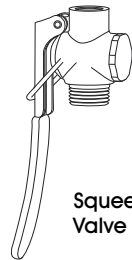Squeeze Valve x 1

## SAFETY

**CAUTION!** In the event of a chemical splash or spill, affected area of the victim's body must be flushed immediately for at least 15 minutes. For chemical spills to the eyes, user should hold eyes open as wide as possible to permit water to reach all areas around the eye.

**CAUTION!** Rinsing affected area for 15 minutes is only the first step in treating exposure to a hazardous chemical. Following rinsing, victim should be examined and treated by a doctor or other trained medical specialist as soon as possible.

## INSTALLATION


**NOTE:** Unit should be installed by licensed plumber.

**NOTE:** Thoroughly clean and flush supply line prior to installing unit. Pipe shavings, scale, tape and other debris can be carried through a pipe and into unit when system is activated. This debris can damage valve components and interfere with proper operation of the valve.

- The water supply line should have outside diameter of at least 1/2" and should be copper or other corrosion-resistant material.

- Unit has a 3/8" NPT male thread on inlet. (See Figure 1) Use Teflon® tape or other pipe sealant to seal thread. Do not apply sealant in a way that will allow it to enter into unit.

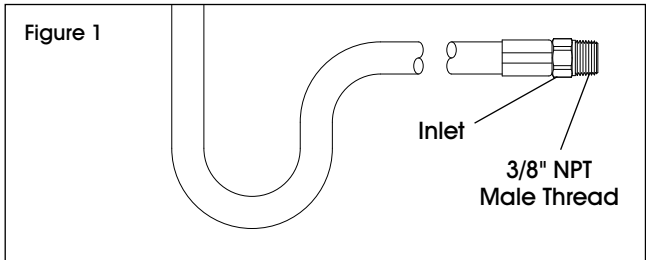## INSTALLATION CONTINUED

**NOTE:** Install unit with adequate clear space around. In an emergency, unit should be readily accessible to the user without interference from other equipment, such as mixing faucets, pegboards, instrumentation, etc.

**NOTE:** Install unit so handle and spray heads face forward toward user.

**NOTE:** No sharp edges, corners or objects should rub against hose. Ensure unit is accessible from wall on which it is installed.

**NOTE:** Unit must pull out easily from wall bracket without restriction. Keep hose free from obstructions to prevent hose from kinking.

## MAINTENANCE

**NOTE:** Unit must be inspected on a regular basis.

**Weekly Activation:** This is required by the provisions of ANSI Z358.1-2014.

**Monthly Maintenance Inspection:** This should incorporate the recommended weekly activation of the unit, as well as a thorough inspection of the unit to ensure that it is in good operating condition and showing no signs of wear. Pay attention to the spray heads, valve and hose assembly.

**Annual Compliance Inspection:** All emergency equipment should be inspected at least annually to verify continued compliance with the provisions of ANSI Z358.1-2014.

Clean unit using only a soft cloth and soapy water. Do not use abrasives, detergents or other cleaners that can damage the finish on the unit.

**WARNING!** Never use cleaning agent that contains ammonia near a laboratory service fitting or emergency eyewash unit, as ammonia will aggressively corrode brass contained in these products.

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