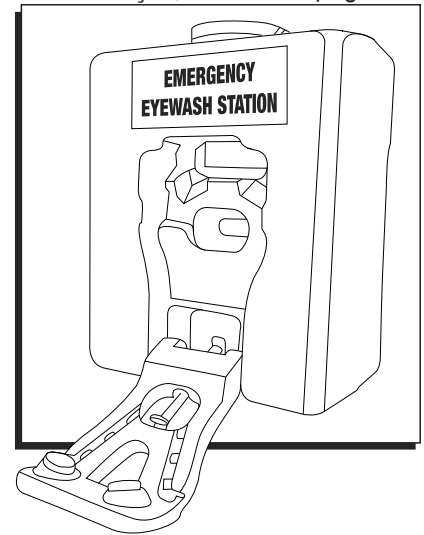
# ULINE H-9832

## SLIM FLOW EYEWASH STATION

1-800-295-5510  
uline.com

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### SAFETY

**WARNING!** Failure to comply with these instructions may result in serious personal injury, serious illness, blindness or death.

- The Slim Flow must be level for proper operation.
- The unit may weigh over 80 lbs. when filled with solution. Assistance may be required when installing unit on the mounting bracket.
- Employees must be instructed in the use of emergency eyewash stations.
- This unit is intended for emergency use by victims of eye contamination until medical treatment is obtained. In all cases, the victim should seek medical attention immediately after flushing.
- An individual or department must be assigned the task of checking all self-contained eyewash units on a weekly basis.
- The use of water preservative is recommended with this eyewash station to prevent the growth of harmful bacteria in the eyewash tank.
- Use of a fluid beyond the expiration date may result in serious personal injury.

### INSTALLATION

#### LOCATION

The ANSI Standard Z358.1-2014 requires that:

1. The Slim Flow should be positioned with the nozzle(s) no less than 33" and no greater than 45" from the floor and 6" minimum from the wall or nearest obstruction.
2. Eyewash units should be in accessible locations that require no more than 10 seconds to reach. The eyewash should be located on the same level as the hazard, and the path of travel should be free of obstructions that may inhibit the immediate use of the equipment. For a strong acid or a strong caustic, the eyewash should be immediately adjacent to the hazard.
3. Each eyewash location should be identified with a highly visible sign positioned so the sign is visible within the area served by the eyewash. The area around the eyewash should be well-lit.
4. The eyewash station must be installed in an easily accessible location in the immediate area of hazard to the employee.

The Slim Flow may be placed on any level surface of required height, such as a table or bench. Allow for proper drainage and disposal of the flushing solution.

The unit may also be mounted on a wall or post using the metal mounting bracket. The unit must be level for proper operation.

## INSTALLATION CONTINUED

### MOUNTING

Using the included hardware, mount bracket to wall stud between 53 and 65" high. Hang eyewash station from the bracket, ensuring spray heads are between 33 and 45" from the floor. (See Figure 1)

**IMPORTANT!** Do not fill eyewash unit before mounting. Filled unit is approximately 80 pounds.

Fill station with clean, potable water to bottom of cap opening, then add the included eyewash preservative and screw on cap. Record the date of filling on the inspection tag.

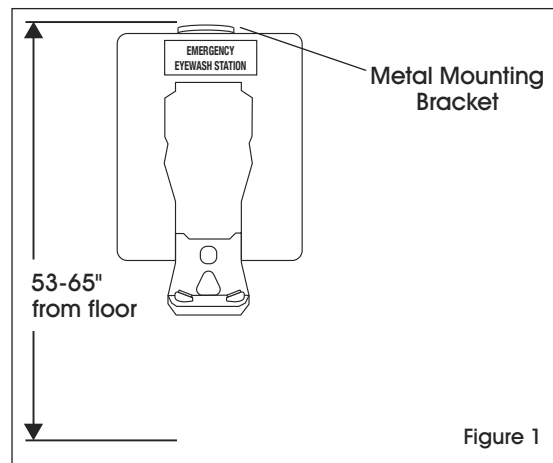
Figure 1

## OPERATION

Employees must be instructed in the use of emergency eyewash stations. Obtaining emergency flushing within the first 10 seconds in eye emergencies is critical. It is required that training be provided to employees in each department where eye hazards exist. This training should be scheduled on a regular basis to keep both current and new employees aware of the location of all eyewash devices, as well as to reinforce the importance of obtaining immediate eye washing in emergencies.

**NOTE:** The sooner the injured person can flush his or her eyes, the less likely there will be of eye damage. Seconds can make the difference in preventing serious eye injury or blindness.

**NOTE:** All employees must be reminded this unit is for emergency use only. Tampering with this unit or using it for non-emergency purposes may cause the unit to not operate properly in an emergency situation.

In an emergency, the eyes may involuntarily clamp shut and the individual may be experiencing pain, anxiety and disorientation. They may need assistance from other workers in getting to and using the eyewash station.

1. Start the fluid flow by pulling down on the black eyewash arm.
2. The injured person should then lower his or her head into the eyewash station toward the nozzles.
3. When eyes are in the stream of fluid, use thumb and forefinger of each hand to open eyelids. Hold eyelids open until thoroughly flushed. 15 minutes is recommended.
4. After flushing, seek medical attention immediately.
5. Safety and supervision personnel must be notified when the unit has been used and must be refilled.

## INSPECTION

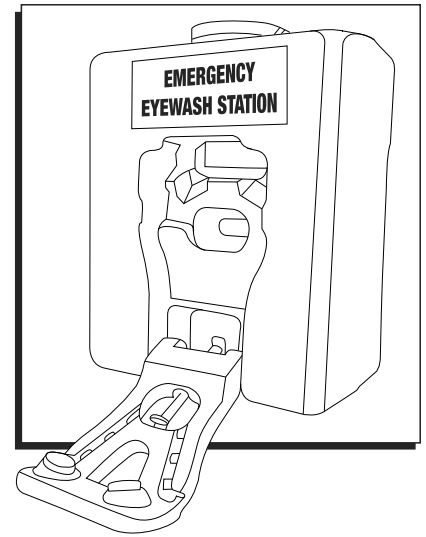
An individual or department must be assigned the task of checking all self-contained eyewash units on a weekly schedule to ensure the unit is in proper operating condition with adequate levels of flushing fluid. Records of these weekly inspections should be kept with the unit. This eyewash station should be cleaned and refilled every six months. Use of a fluid beyond the expiration date may result in serious personal injury.

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