
ULINE H-4737, H-4738
H-4739
OHAUS EXPLORER®
BALANCE SCALE


1-800-295-5510
uline.com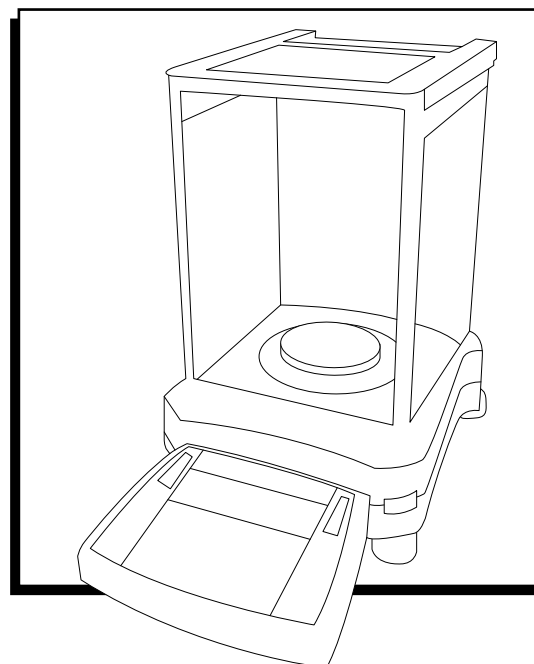

TROUBLESHOOTING

OPERATING ISSUES	POSSIBLE CAUSES	RECOMMENDATIONS
Balance will not turn on.	No power to balance.	Verify connection and voltage.
Poor accuracy.	Improper calibration. Unstable environment.	Perform calibration. Move balance to suitable location.
Cannot calibrate.	Calibration menu is locked. Unstable environment. Incorrect calibration masses.	Turn calibration menu lock off. Move balance to suitable location. Use correct calibration masses.
Cannot change menu settings.	Sub-menu is locked.	Unlock sub-menu.
Low reference weight.	Reference weight too small. The weight on the pan is too small to define a valid reference weight.	Increase sample size.
Invalid piece weight.	Average piece weight is too small.	Increase average piece weight.
Operation time out.	Weight reading is not stable.	Move balance to suitable location.
Screen displays "-----".	Busy (tare, zero, printing).	Wait until completion.
Scale will not weigh or is displaying inaccurate weights.	Scale has fallen out of calibration.	Recalibrate scale. Refer to page 5 of instructions .

If the troubleshooting section does not resolve your problem, contact Uline Customer Service at 1-800-295-5510.

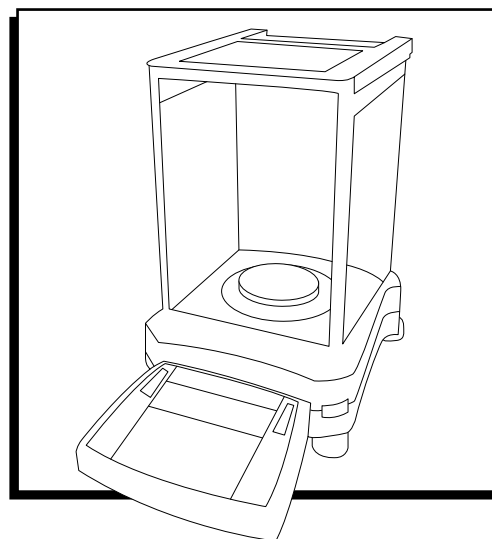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