
ULINE H-5276, H-5278
H-5279, H-6097
OHAUS ADVENTURER®
BALANCE SCALE


1-800-295-5510
uline.com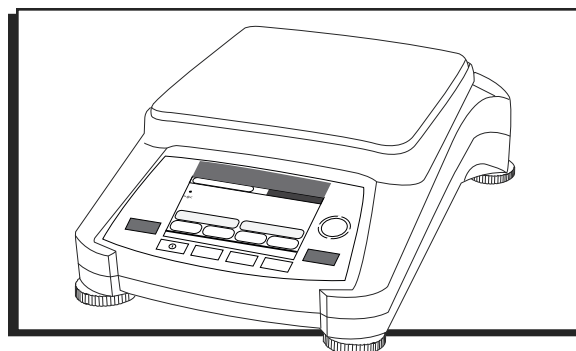

TROUBLESHOOTING

OPERATING ISSUES	POSSIBLE CAUSES	RECOMMENDATIONS
Balance will not turn on.	No power to balance.	Verify connection and voltage.
Poor accuracy.	Improper calibration. Unstable environment.	Platform calibration. Move balance to suitable location.
Cannot calibrate.	Calibration menu locked. Approved Mode set to "on". Unstable environment. Incorrect calibration masses.	Turn Calibration menu lock off. Turn Approved Mode off. Move balance to suitable location. Use correct calibration masses.
Cannot change menu settings.	Sub-menu locked. Approved Mode set to "on".	Unlock sub-menu. Turn Approved Mode off.
Low Reference weight.	Reference weight too small. The weight on the pan is too small to define a valid reference weight.	Increase sample size.
Invalid Piece Weight.	Average piece weight is too small.	Increase average piece weight.
Operation timeout.	Weight reading is not stable.	Move balance to suitable location.
Screen displays "-----".	Busy (tare, zero, printing, waiting for a stable weight).	Wait until completion.
Scale will not weigh or is displaying inaccurate weights.	Scale has fallen out of calibration.	Recalibrate scale. Refer to page 3 of instructions .

If the troubleshooting section does not resolve your problem, contact Uline Customer Service at 1-800-295-5510.

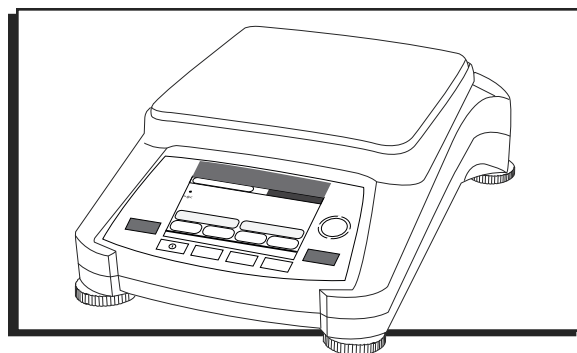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