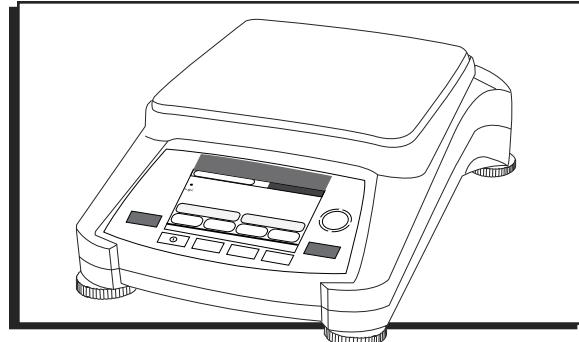
ULINE H-5276, H-5278
H-5279, H-6097

OHAUS ADVENTURER®
BALANCE SCALE

1-800-295-5510
uline.com


TROUBLESHOOTING

OPERATING ISSUE	CAUSES	RECOMMENDATIONS
Balance will not turn on.	No power to balance.	Verify connection and voltage.
Poor accuracy.	Improper calibration. Unstable environment.	Platform calibration. Move balance to suitable location.
Cannot calibrate.	Calibration menu locked. Approved Mode set to "on". Unstable environment. Incorrect calibration masses.	Turn Calibration menu lock off. Turn Approved Mode off. Move balance to suitable location. Use correct calibration masses.
Cannot change menu settings.	Sub-menu locked. Approved Mode set to "on".	Unlock sub-menu. Turn Approved Mode off.
Low Reference weight.	Reference weight too small. The weight on the pan is too small to define a valid reference weight.	Increase sample size.
Invalid Piece Weight.	Average piece weight is too small.	Increase average piece weight.
Operation timeout.	Weight reading is not stable.	Move balance to suitable location.
Screen displays "-----".	Busy (tare, zero, printing, waiting for a stable weight).	Wait until completion.

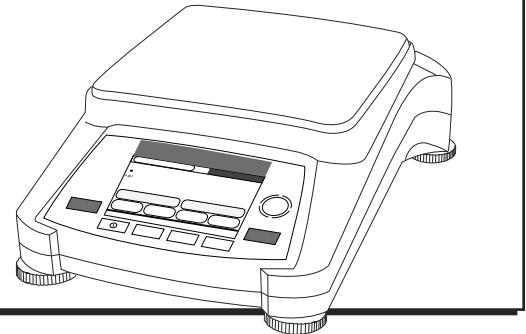If the troubleshooting section does not resolve your problem, contact Uline Customer Service at 1-800-295-5510.

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