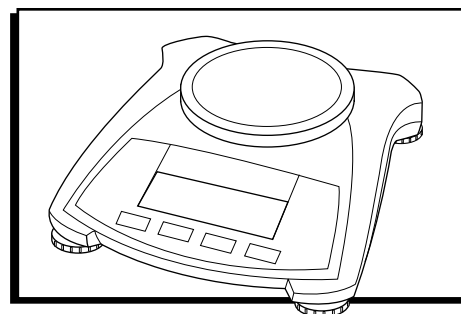
ULINE

**OHAUS SCOUT®
BALANCE SCALE**

1-800-295-5510
uline.com

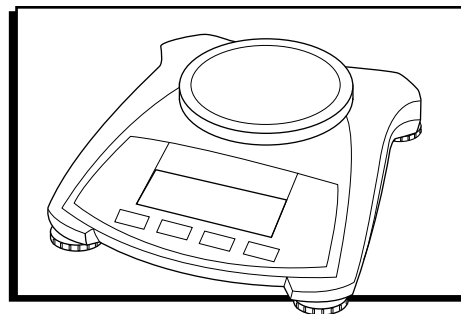
TROUBLESHOOTING

OPERATING ISSUES	RECOMMENDATIONS
Cannot turn on.	No power to balance. Bad power source. AC adapter or batteries. Faulty keypad.
Poor accuracy.	Improper calibration; Unstable environment.
Cannot calibrate.	Security switch in locked position.
Scale will not weigh or is displaying inaccurate weights.	Move transportation lock to unlocked position. Recalibrate scale.

ERROR CODE	RECOMMENDATIONS
<i>Err 8.1</i>	Weight exceeds power on zero range. Remove weight, restart and recalibrate.
<i>Err 8.2</i>	Weight below power on zero range. Platform may have been removed before turning on the scale. Replace platform, start the scale and recalibrate.
<i>Err 8.3</i>	Overload (weight exceeds rated capacity).
<i>Err 8.4</i>	Under load (platform removed).
<i>Err 8.5</i>	Tare out of range.
<i>Err 8.6</i>	Displayed value exceeds 999999 (possible in totalization mode).
<i>rEF.Err</i>	Parts counting or percentage error – sample weight <1d. Balance shows error then exits parts counting or goes to <i>Clr.APU</i> .
<i>Lo.rEF</i>	Percent Reference weight or APW is too low for accurate results.
<i>CLL E</i>	Fail to do calibration.
<i>USB.Err</i>	Cannot find menu or application file in U-disk.

If the troubleshooting section does not resolve your problem, contact Uline Customer Service at 1-800-295-5510.

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