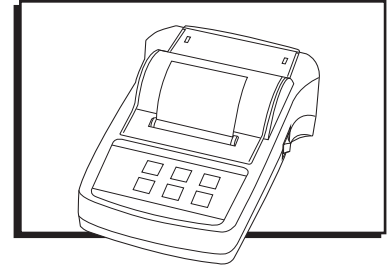
ULINE H-8093
OHAUS® SF40A
SCALE PRINTER

1-800-295-5510
 uline.com

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TROUBLESHOOTING

OPERATING ISSUE	CAUSES	RECOMMENDATIONS
Printout unreadable.	Ribbon worn or tangled. Service life of printing unit has ended.	Change ribbon cartridge or tense the ribbon properly. Contact Uline Customer Service at 1-800-295-5510.
Pilot lamp does not light.	No power. Power supply fault.	Switch printer on. Contact Uline Customer Service at 1-800-295-5510.
Pilot lamp flashes.	No connection printer to device. Printer unable to recognize device settings for the connected scale.	Make sure that the printer and the scale are connected. Switch the printer off/on. Alternatively change settings of the printer or the connected scale. (See page 4 of instructions)
----- Timeout -----	Data transfer not completed within 45 seconds.	Move scale to a more stable location. Check the scale.
----- Over Load -----	Balance in overload range. Zeroing out of the zero range.	Check the sample weight. Change the zero range of the balance. Change the setting of the printer from zero to tare.
----- Under Load -----	Scale in underload range. Weighing pan is not in position.	Check the scale.
---- Transmission Error----	Scale does not recognize printer instruction.	Check the interface parameter settings on the printer and scale.
----- Not Executable -----	Scale cannot execute printer instruction.	Check the scale.

TROUBLESHOOTING

OPERATING ISSUE	CAUSES	RECOMMENDATIONS
--- Date/Time Not Set --- ----- Memory Error -----	Printer error messages.	Check the scale. Reset the printer. Set date/time. Backup battery depleted. Contact Uline Customer Service at 1-800-295-5510.
----- Memory Full -----	Statistics or Totaling Memory is full.	Start a new Statistics or Totaling application.
-- Load/Unload Weight --	No sample or the same sample is on the weighing pan.	Put a new sample on the weighing pan or remove it and put sample on the weighing pan again.
Unit change not allowed!	Changed unit during Statistics or Totaling application.	Reset the scale. Change the unit back to the original unit setting.
----- Out of range -----	Sample weight is out of range.	Sample weight must be in the range of 70-130% of the current average value.

If the troubleshooting section does not solve your problem, contact Uline Customer Service at 1-800-295-5510.

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