
ULINE H-9884, H-9885
 H-9886
BALANCE SCALE


1-800-295-5510
 uline.com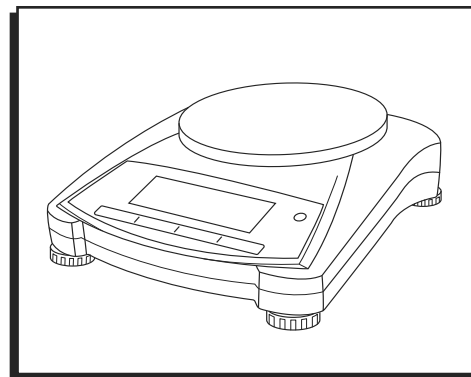

TROUBLESHOOTING

OPERATING ISSUE	POSSIBLE CAUSE	REMEDY
Cannot turn on	No power to scale.	Verify connections and voltage.
Poor accuracy	Unstable environment. Improper calibration.	Move scale to suitable location away from airflow and vibrations. Perform calibration.
Cannot calibrate	Unstable environment. Incorrect calibration weight.	Move scale to suitable location away from airflow and vibrations. Use correct calibration weight.
Cannot access mode	Mode not enabled.	Enter menu and enable mode.
Cannot access unit	Unit not enabled.	Enter menu and enable unit.
<i>Lo rEF</i>	Reference weight is too low.	Increase reference weight.
<i>rEF Err</i>	Parts counting – sample weight < 1d.	Shows error, exits mode or goes to <i>Clr.RPU</i> .
<i>Err 3.0 CAL</i>	Incorrect calibration weight.	See Calibration on page 3.
<i>ERR 4.4 FULL</i>	RS232 Buffer is full.	Set Handshake on, see Interface User Manual.
<i>Err 8.1 LoRd</i>	Power on zero range exceeded.	Clear pan; check shipping lock setting.
<i>Err 8.2 LoRd</i>	Power on zero under range.	Install pan; check shipping lock setting.
<i>Err 8.3 LoRd</i>	Overload.	Load exceeds scale maximum capacity.
<i>Err 8.4 LoRd</i>	Under load.	Reading below min. range; re-install pan.
<i>Err 8.6 999999</i>	Displayed value > 999999.	Result exceeds display capability.
<i>Err 9 dRtR</i>	Internal data error.	Contact Uline Customer Service.
<i>Err 13 MEM</i>	Fail to write EEPROM.	Contact Uline Customer Service.
<i>Err 53 CSUM</i>	Invalid checksum data.	Contact Uline Customer Service.

If troubleshooting section does not resolve or describe problem, contact Uline Customer Service at 1-800-295-5510.

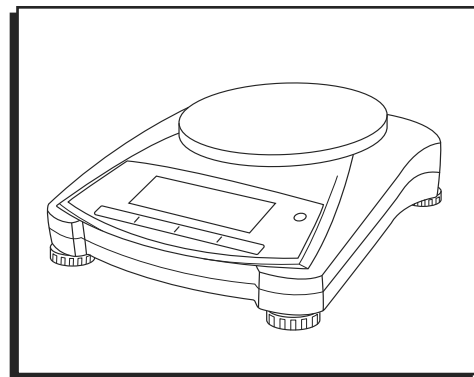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